


抗体薬物複合体 (ブレンツキシマブ ベドチン)
Antibody-Drug Conjugates (Brentuximab vedotin)

C240226B

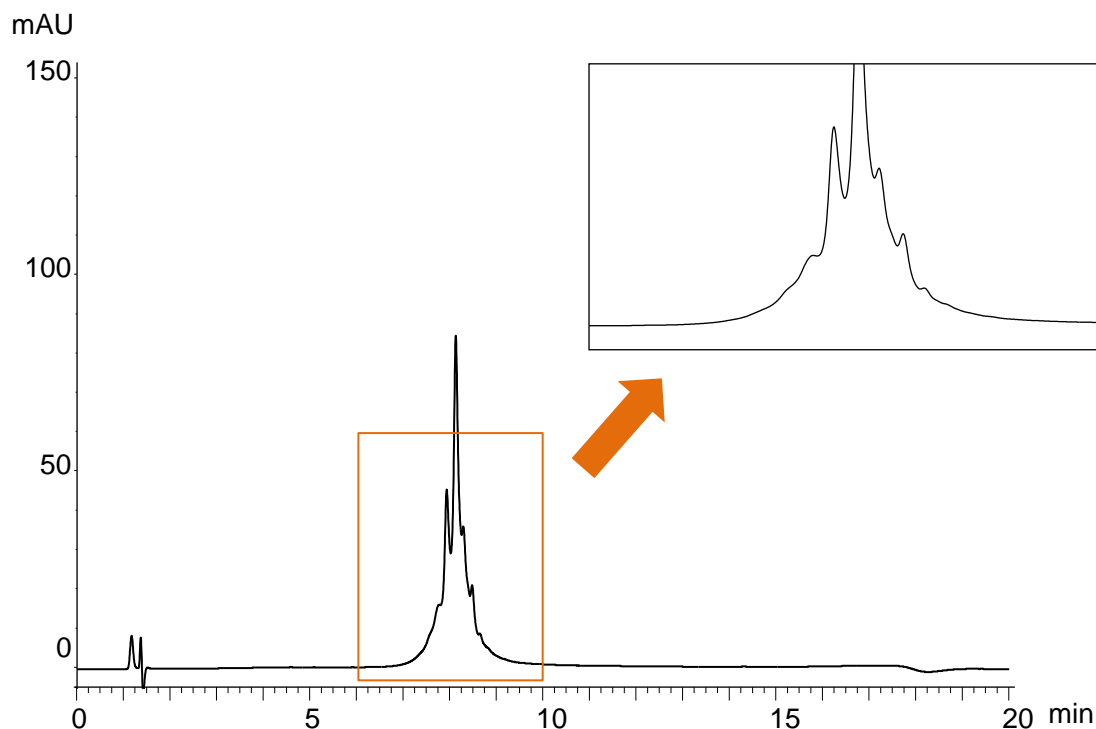

Column : Accura BioPro IEX SF (3 μ m)
100 X 4.6 mm I.D.

Eluent : A) 10 mM MES*-NaOH (pH 6.0)
B) 10 mM MES*-NaOH (pH 6.0) containing 1.0 M NaCl
0–20%B (0–15 min)

Flow rate : 0.5 mL/min

Temperature : 25°C

Detection : UV at 280 nm


Injection : 10 μ L (1 mg/mL)

Sample : Brentuximab vedotin for injection
Courtesy of Prof. S. Manabe, Hoshi University, Japan/Tohoku University, Japan

*2-(*N*-morpholino) ethanesulfonic acid

抗体薬物複合体 (ブレンツキシマブ ベドチン)
Antibody-Drug Conjugates (Brentuximab vedotin)

AK240328C

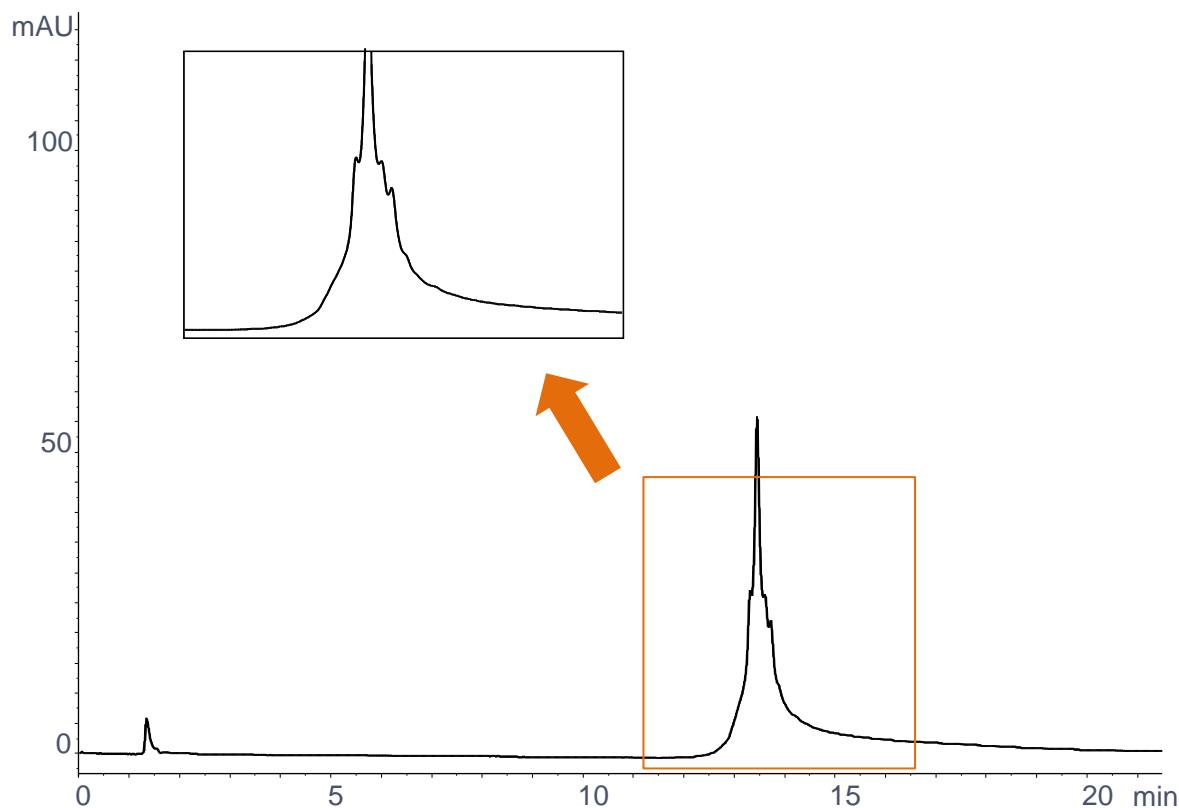

Column : Accura BioPro IEX SF (3 μ m)
100 X 2.1 mm I.D.

Eluent : A) 20 mM CH₃COONH₄-CH₃COOH (pH 5.6)
B) 140 mM CH₃COONH₄-10 mM NH₄HCO₃ (pH 7.4)
0%B (0–2 min), 0–100%B (2–18 min), 100%B (18–22 min)

Flow rate : 0.1 mL/min

Temperature : 25°C

Detection : UV at 280 nm

Injection : 2 μ L (1 mg/mL)

Sample : Brentuximab vedotin for injection

Courtesy of Prof. S. Manabe, Hoshi University, Japan/Tohoku University, Japan